

Technical Bulletin





Author: Larry Pastwik - Product Specialist, Analytical Instruments

Date: 30 November 2004

MARK II CONTACT CLEANING

Following is the procedure for cleaning the electrical contact points within the ABBE MARK II PLUS Refractometer:

1. Unplug instrument from power source.
2. Remove large shadow line adjusting knob by loosening set screw located through cutout in knob edge.
3. Remove right side cover after removing screws located at the bottom of the panel.
4. Clean contact surfaces with a cotton swab and Isopropyl Alcohol, Methanol or window cleaning solution.
5. Reassemble, plug in and test.
6. Repeat process if required.

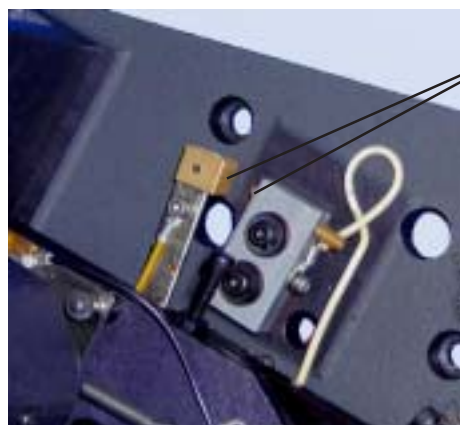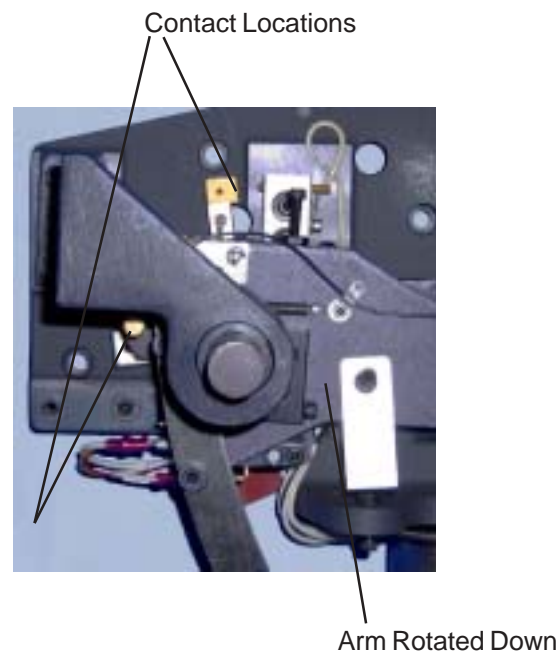

Contact Surfaces

